

	BIO-PM 4.30 Rainin EDP and Pipet-lite Pipettors	
	<i>Document #: 1938</i>	<i>Page 1 of 1</i>
	<i>Revision #: 1</i>	<i>Issued Date: 11/14/2017</i>
	<i>Document Manager: Kristin Schelling</i>	<i>Approved By: Jeffrey Nye</i>

4.30 Rainin EDP And Pipet-lite Pipettors

4.30.1 Scope

To provide an accurate measurement of very small quantities with a minimum of repetitive motion trauma to the operator.

4.30.2 Background Information

Owner's manuals and reference guides for Pipet-Lite and EDP3.

4.30.3 Maintenance

4.30.3.1

Read all instructions completely before use.

4.30.3.2

The Pipet-Lite has an autoclavable shaft and tip-ejector. No other parts are autoclavable.

4.30.3.3

These pipettors must be sent in for annual calibration to maintain traceable laboratory standards.

Decontaminate these instruments in the following manner before shipping to Rainin:

4.30.3.3.1

Electronic models: Wipe down the outside and the tip mounts with 10% bleach solution and allow to air dry.

4.30.3.3.2

Manual models: In addition to autoclaving the shaft and tip-ejector, other parts may be wiped with ethanol, then dried. See owner's manual.