

	BIO-PM 4.31 Decontamination of Regular Rainin Pipetman Pipettors Prior to Annual Recalibration	
	Document #: 1939	Page 1 of 1
	Revision #: 1	Issued Date: 11/14/2017
	Document Manager: Kristin Schelling	Approved By: Jeffrey Nye

4.31 Decontamination Of Regular Rainin Pipetman Pipettors Prior To Annual Recalibration

4.31.1

Remove tip ejector.

4.31.2

Unscrew blue ring around shaft. Remove.

4.31.3

Carefully remove shaft. Try to keep O-rings and seals on piston and piston assembly together.

4.31.4

Soak all parts in a 10% bleach solution for about 15 minutes.

4.31.5

Thoroughly rinse all pieces with deionized water. Shake off excess water.

4.31.6

Rinse all parts with 70% ethanol. Shake off excess.

4.31.7

Place parts in a heat-resistant container and put the whole thing in the vacuum drying oven set on 1.

4.31.8

Keep in the oven till dry. Let cool and reassemble.

4.31.9

Package and ship UPS 2 day to Rainin. Insure each pipettor for \$250.00.