

	BIO-PM 4.5 Ohaus Balance	
	<i>Document #: 1948</i>	<i>Page 1 of 1</i>
	<i>Revision #: 1</i>	<i>Issued Date: 11/14/2017</i>
	<i>Document Manager: Kristin Schelling</i>	<i>Approved By: Jeffrey Nye</i>

4.5 Ohaus Balance

4.5.1 Scope

To ensure that the balance is operating correctly. The balance shall be calibrated annually, or at the time of any service or repair by a certified technician. The calibration performed by the certified technician shall serve as the performance check.

4.5.2 Background Information

See operator's manual

4.5.3 Operation

4.5.3.1

Press the on/off button to turn the balance on. The display will subsequently indicate zero (weighing mode).

4.5.3.2

Load item on the pan to be weighed. Wait for the balance to stabilize and record reading.

4.5.3.3

For taring, place container on the pan, press the re-zero button to initiate the tare cycle. When the display indicates zero the container can be used to weigh the amount of sample.

4.5.3.4

To turn the balance off, press the on/off button. The balance should be left on unless being serviced or the balance is being moved.